


User Manual

Computer Program „BOECO WINPC”

Manual number:
01 2015-07-ENG

Version: 5.X.X

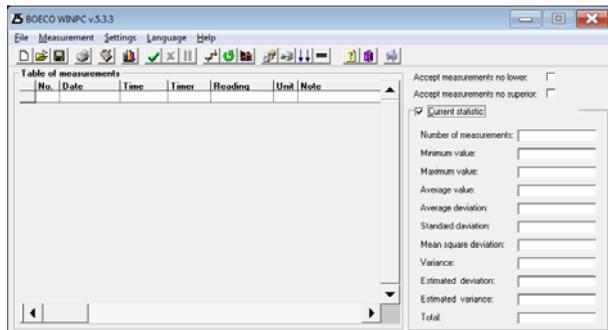

BOECO
Germany

Boeckel + Co (GmbH + Co) KG
Roedingsmarkt 33
20459 Hamburg
Germany
Phone: + 49 40 325627-0
Fax: + 49 40 325627-41
E-mail: info@boeco.com
Internet: www.boeco.com

TABLE OF CONTENTS

1. INTENDED USE	3
2. INSTALLATION	3
2.1. System requirements.....	3
2.2. Installation procedure	4
2.2.1. Program update	6
3. STARTING PROGRAM	6
3.1. Main window.....	6
3.1.1. Window description.....	7
3.1.2. Close the program	9
3.1.3. Help.....	10
4. SETTINGS MENU	10
4.1. Options	10
4.2. Cycle.....	13
4.3. Chart.....	14
5. FILE MENU	15
6. MEASUREMENTS	17
6.1. Automatic operation.....	17
6.2. Read	17
6.3. Tarring	18
6.4. Continuous transmission	18
6.5. Reading non-stable data	19
6.6. Calibration report	20
6.7. Chart.....	21
7. LANGUAGE	22

1. INTENDED USE

Program „ **BOECO WINPC**” has been designed for collecting measurement data from scales made by **BOECO** and performing some statistical calculations.

Basic program functions:

- Reading data from scales, recording measurements in a file,
- Possibility of manual or automatic readings of measurements,
- archiving of measurement data in a file,
- visualization of measurement data in a chart,
- statistical processing of measurement data,
- creating reports from the selected set of measurements.

Notice:

Boeco shall not be liable for any errors resulting from wrong use of the program.

2. INSTALLATION

2.1. System requirements

Minimum requirements:

- Processor INTEL or AMD \geq 1200 MHz,
- RAM minimum 256 MB (recommended 512 MB),
- Graphic card SVGA (resolution minimum 1024x768 with 24-bit colour palette),
- Communication port RS-232,
- Free space on a Hard Drive \geq 1 GB,
- Mouse,
- Keyboard,
- A colour monitor with a resolution of 1024x768,
- System Windows XP/2000/Vista/7,
- Printer (format A4),
- Software for reading pdf files (for manuals),
- Basic computer operation skills.

2.2. Installation procedure

In order to install the program put the installation CD to the CD-ROM drive. An installation program of „**BOECO WINPC**” should start running automatically. If the installation is to be performed manually run program **<Install_BOECO_WINPC_5.x.x.exe>** in the main folder of the CD.


Follow the instructions in the subsequent windows.

The welcome window of the installation wizard:




Press „**Next**” to continue.

The Ready to Install window:

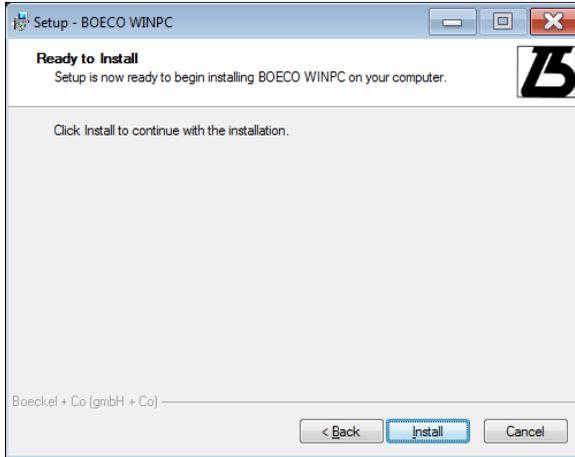

Press button „**Install**” to continue.

The installation completion window:

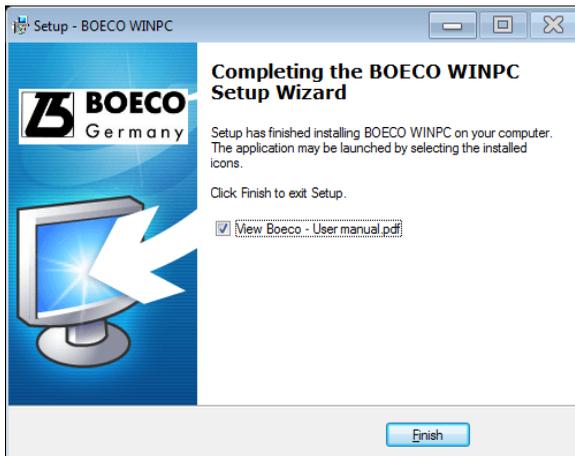

Then „**BOECO WINPC**” will be quickly installed.
Press „**Finish**” to close the installation wizard.

2.2.1. Program update

In case when program „ **BOECO WINPC**” has already been installed on a computer and when the user starts to install a new version of the program, the whole installation process is similar to that when it is installed for the first time.

Notice:

Before the installation wizard is started it is advisable to make a backup copy for the program folder. Default folder is:

<C:\Program Files (x86)\Boeco\BOECO WINPC>.

3. STARTING PROGRAM


The program may be run by clicking the shortcut icon on the desktop:

The program may also be started from the START menu: „**START/Programms/Boeco/BOECO WINPC**”.



3.1. Main window

After starting the program the main window is displayed:

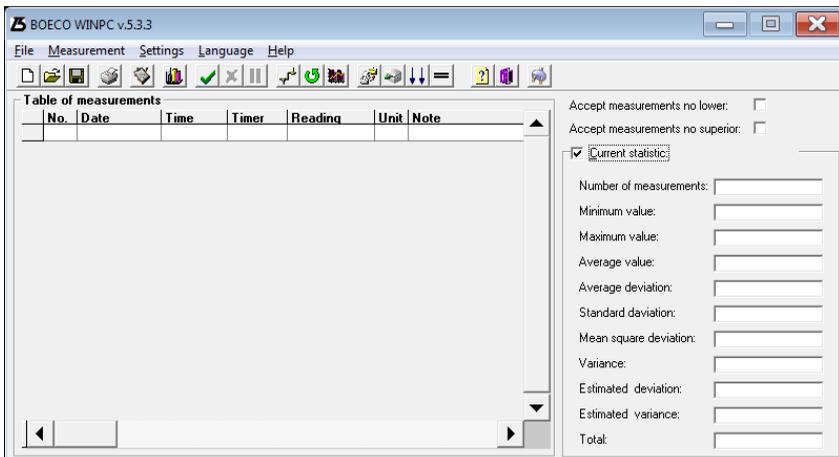

It allows for displaying collected weighings and statistics in real time.

The main program options can be accessed from the menu or buttons in the tool bar. There are additional buttons to set data to be accepted by the program:

- If the cyclical mode is enabled the icon  is displayed, otherwise  can be seen, the cycle interval is set in the program which requests a reading in every cycle.
- If the continuous mode is enabled the icon  is displayed, otherwise  can be seen,
- If the scales sends stable data it is signalled by icon , otherwise icon  can be seen.

Notice:


*When the weighings are collected the statistics are calculated in real time after every weighing. It can cause some problems on computers with low processor capacity. The program may slow down some operations or even the operation in real time may be impossible. In such case option „**Current statistics**” needs to be disabled.*

3.1.1. Window description

Title bar – displays program name and version. Standard windows buttons allows for minimizing, maximizing or closing the window:



Main menu – a menu with all functions of the program:

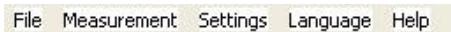

Buttons in the toolbar – buttons with direct access to some functions:



Buttons' description:

	Open new series of measurements
	Open an existing series of measurements in file with extension *.ser
	Saves current weighings from the table to a file
	Prints current measurements and statistics
	Opens the report window
	Shows the chart of weighings
	Starts automatic operation
	Terminates automatic operation
	Holds automatic operation
	Opens program settings
	Opens cycle settings
	Opens chart settings
	Reads one measurement from a scale
	Tares scale
	Initiates or terminates continuous transmission
	Enables / disables non-stable readings
	Displays help
	Closes program

Table of measurements – an area for displaying measurements:

Table of measurements						
No.	Date	Time	Timer	Reading	Unit	Note
1	2010-09-13	15:00:55	0	0,990	kg	
▶ 2	2010-09-13	15:00:55	0	0,990	kg	

Window of Statistics – displays statistical information about measurements.

Accept measurements no lower:

Accept measurements no superior:

Current statistic:

Number of measurements:

Minimum value:

Maximum value:

Average value:

Average deviation:

Standard deviation:

Mean square deviation:

Variance:

Estimated deviation:


Estimated variance:

Total:

After choosing one of options „**Accept measurements no lower / no superior**”, a window to describe a limit value will appear.

3.1.2. Close the program

Closing the application is possible after pressing button  or choosing item „**File/Exit**” in the menu:

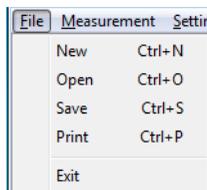

Closing the program and some windows is preceded by displaying a dialogue window to confirm or contradict the operation.:



3.1.3. Help


Info window will be displayed after pressing  or choosing item „**Help**” in the menu:




4. SETTINGS MENU

4.1. Options

The access to menu item „**Options**” is possible after pressing button  or entering submenu „**Settings/Options**”:

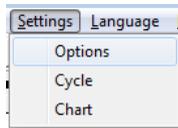

In menu „**Options**” parameters of communication with a scale are set. Connect the scale to an RS232 COM port or Ethernet and then carry out configuration of parameters according to the description.

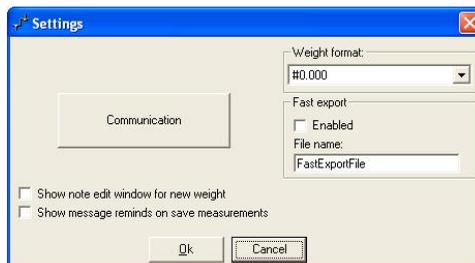

Communication

The button for accessing a configuration window for defining the connection with a scale.


- **Communication via RS-232**




Description:

Port Number Serial port number to be used for communication with scales

Baud rate Communication speed, the same as set in scales

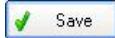


Save settings

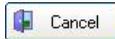

Cancel settings


- **TCP communication**

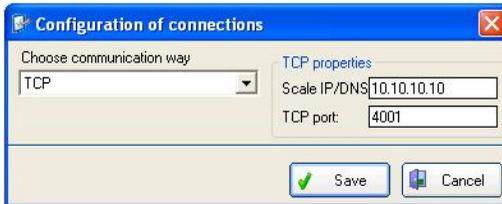

Description:

Scale IP/DNS IP address used by a scale

TCP port TCP port number used by a scale

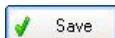


Save settings



Cancel settings

Measurement Format

Set displaying format.

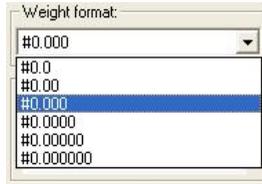

After modifying it the changes will be visible after restarting the program in the table of measurements.

Fast export


After enabling the option measurements from the table will be automatically saved to file *.csv with given name. Subsequent savings are added to the file.



Choosing the item „**Save**” in menu „**File**” causes clearing the table and saving measurements in the right file. Saving measurements to a file with extension *.ser is not possible then.

Editing description

When option **<Show note edit window for new weight>** is selected, after every measurement a window can be displayed for typing a description of the measurement.

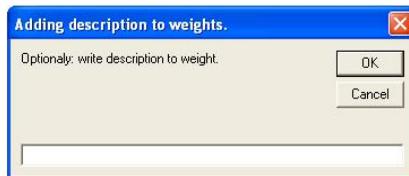


Editing records from the table of weighings is possible as well (main program window) after double-click on a measurement.

Saving remainder


When option **<Show message reminds on save measurements>** is selected then when 18 measurements are overpassed, so the list in the table of measurement is full, the message window will be displayed. All changes in configuration will be introduced after pressing button „OK”.

4.2. Cycle

Access to option „**Cycle**” is possible after pressing button  or entering the menu item „**Settings/ Cycle**”:

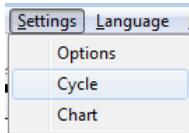

In the program you can set a time interval between subsequent reading and enable the cyclical readouts.

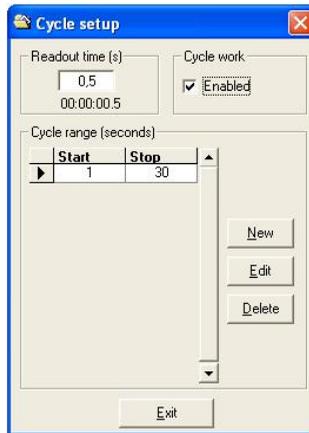

Subsequent readouts with time interval (s)


Setting the time in seconds for subsequent readings of measurement from a scale. The minimum range is 0.5 s – and maximum 86399 s.

Cyclical operation


Enables operation of the program with the declared step. Numbers in fields mean quantities of seconds which have elapsed from the first measurement.

4.3. Chart

Setting options for „**Chart**” is accessible after pressing  or choosing menu item „**Settings/Chart**”:

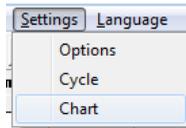


Window „**Chart setup**” allows for changing the look of the chart:

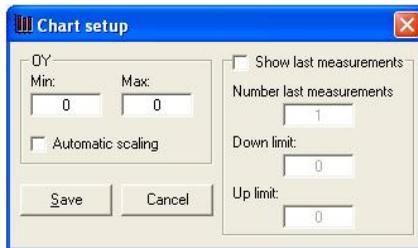

In fields **Min** and **Max** set the range of values you want to see in the chart. If you do not want to limit the range tick the option „**Automatic scaling**”.

After enabling option „**Show last measurements**”, you can limit the number of measurements put in the chart to the quantity declared in field „**Number last measurements**”. Additionally you can display two horizontal lines that would represent some limitations for measurements or simplify readings.

Example chart:

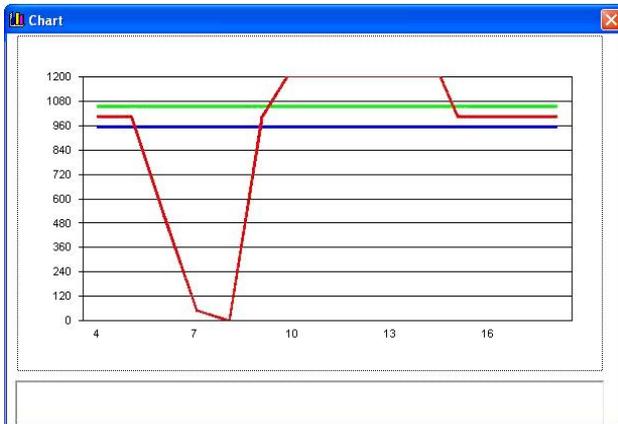

Colours:


- Red** Measurements
- Green** Top threshold
- Blue** Bottom threshold

5. FILE MENU

Different file functions are accessible after pressing buttons:




or choosing items in menu „File”:

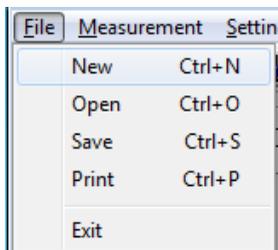

File menu items:

- New** Starting a new series of measurements, deleting current measurements and statistics.
- Open** After opening a file measurements from the file will be recovered and as well as statistics saved in file with extension *.ser. If current measurements are included in the table of measurements a warning is displayed and users can save them before erasing.
- Save** Saving current measurements from the table of measurements. Type a file name and choose a folder to save the file. The saved files will have extensions *.ser (series) or *.csv, depending on settings. After savings the table of measurements is emptied and statistics zeroed.
- Print** Printing weighings and statistics. After choosing the item a printout view is displayed. Press the printer icon in the window to print the data.
- Exit** Closing the program

The view of general report generated from measurements and statistics:

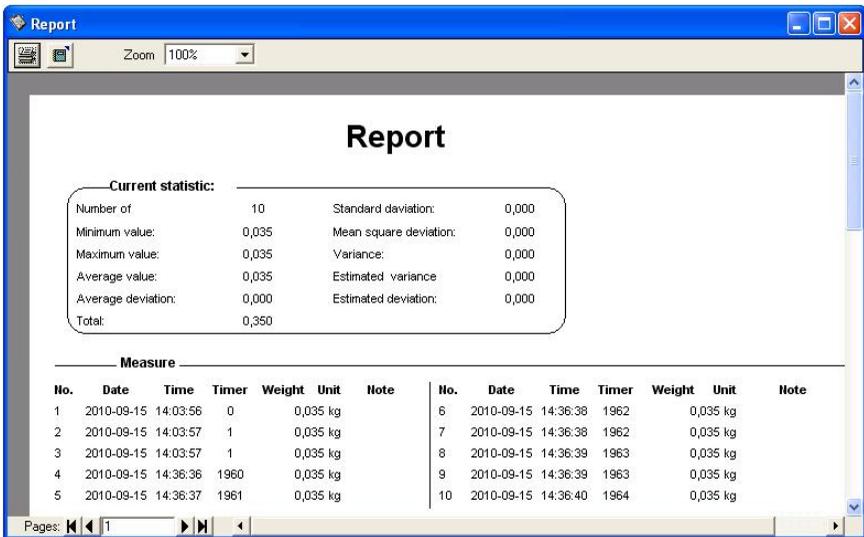

The screenshot shows a software window titled "Report" with a blue title bar and standard Windows window controls. The window contains a report with the following content:

Report

Current statistic:

Number of	10	Standard deviation:	0,000
Minimum value:	0,035	Mean square deviation:	0,000
Maximum value:	0,035	Variance:	0,000
Average value:	0,035	Estimated variance	0,000
Average deviation:	0,000	Estimated deviation:	0,000
Total:	0,350		

Measure

No.	Date	Time	Timer	Weight	Unit	Note	No.	Date	Time	Timer	Weight	Unit	Note
1	2010-09-15	14:03:56	0	0,035	kg		6	2010-09-15	14:36:38	1962	0,035	kg	
2	2010-09-15	14:03:57	1	0,035	kg		7	2010-09-15	14:36:38	1962	0,035	kg	
3	2010-09-15	14:03:57	1	0,035	kg		8	2010-09-15	14:36:39	1963	0,035	kg	
4	2010-09-15	14:36:36	1960	0,035	kg		9	2010-09-15	14:36:39	1963	0,035	kg	
5	2010-09-15	14:36:37	1961	0,035	kg		10	2010-09-15	14:36:40	1964	0,035	kg	

Pages: 1


6. MEASUREMENTS

Notice:

In case of lack of communication between the program and a scale check the configuration of connection in the menu: „Settings/Options/Communication”.

6.1. Automatic operation

The automatic operation is started by pressing button  or choosing item „**Start**” in the menu „**Measurements**”:

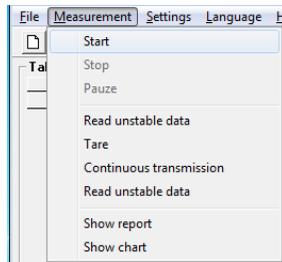


The frequency of readings is set in window „**Cycle setup**” (see section 4.2). If the cyclical operation is enabled the program fetches data in the predefined cycles. In other cases measurements are fetched continuously (with maximum speed for the baud rate). The automatic operation stops after pressing button  or item „**Stop**” in the menu “**Measurements**”.

The automatic operation can be terminated after pressing button  or item „**Pauze**” in the menu “**Measurements**”.

6.2. Read


In order to read data from a scale press button  or choose item „**Read stable data**” in the menu „**Measurements**”:

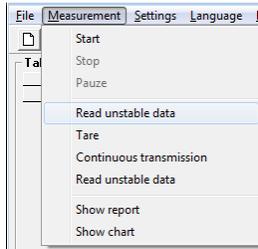

You can press button „Print” on the scale.

6.3. Tarring

A scale cooperating with “Boeco” software can be tarred by pressing button  or choosing item „Tare” in the menu „Measurements”:

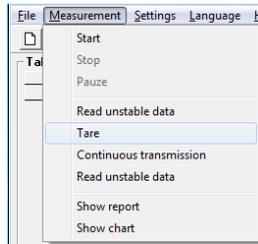


The scale will display **0**, and all performed measurements will be of net value.

6.4. Continuous transmission

Starting the continuous transmission causes the scale to send data to the computer without any requests.

Press button  to start the continuous transmission or choose item „Continuous transmission” in the menu „Measurements”:

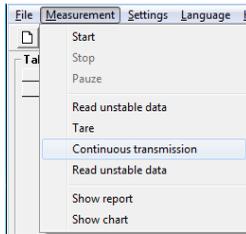


A window informing about receiving data is displayed.



Press button  to stop continuous transmission or press „**Esc**” on the keyboard. Then measurements will be copied to a database.

6.5. Reading non-stable data

If we want to register data without checking if results are stable or not press button  or choose item „**Read unstable data**” in menu „**Measurements**”:

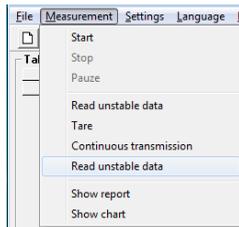


If a scale sends „**non-stable**” data icon  is displayed in the toolbar.


If the scale sends „**stable**” data icon  is displayed in the toolbar.

6.6. Calibration report

In order to show all data send by a scale (not only measurements) press icon  or choose item „**Show report**” in the menu „**Measurements**”. It can be used e.g. for receiving and saving calibration reports.

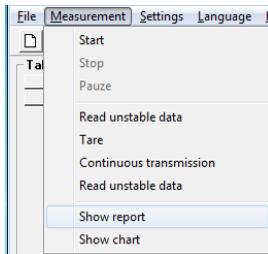

Calibration report window:

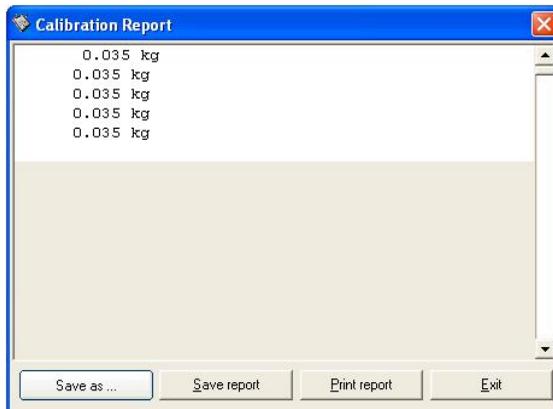

Data can be saved to a text file by pressing button „**Save as...**” and/or print to a *.pdf file by pressing button „**Print report**”.

6.7. Chart

Collected measurements can be presented as a chart. Press button or „**Measurements**” in the menu and submenu „**Show chart**” to draw the chart:


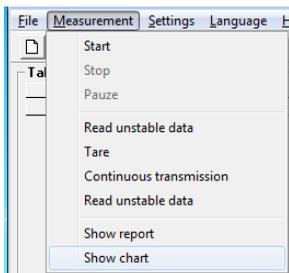


Chart parameters can be changed in window „**Chart setup**” (see section 4.3). In order to print the chart right-click the chart and choose „**Print**” from the pop-up menu.



7. LANGUAGE

Users can change the language after choosing „**Language**” in the menu:

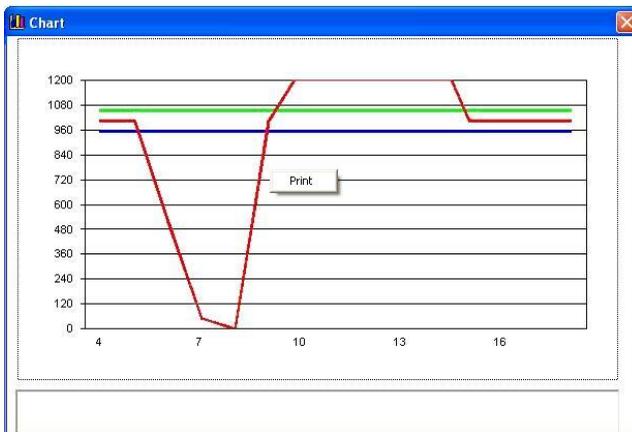

Confirm the choosen language by pressing „**OK**”, then the following message window appears:



After pressing „**OK**” the program needs to be rerun to use the new settings.



Boeckel + Co (GmbH + Co) KG

Roedingsmarkt 33

20459 Hamburg

Germany

Phone: + 49 40 325627-0

Fax: + 49 40 325627-41

E-mail: info@boeco.com

Internet: www.boeco.com